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Insert USB Flash Drive
into USB slot on the
MAC.

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Insert USB Flash Drive
into USB slot on the
PC

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Quick Start Guide

Before you Begin:

Your Moxi Z comes with a partial battery charge and can be used immediately. Alternatively, you can charge the Moxi by connecting it to a computer or power adapter via the supplied USB cable. Be sure to disconnect the USB cable prior to running samples.

Required Materials:

- Moxi Z cassettes and diluent.
- Biological cell sample.
- Pipette and tips for 75µL aliquot.

Sample Prep Options:

-Type S Cassettes. Size range: 3 - 20µm avg. dia. Concentration range: 3,000 to 2,500,000 cells/ml
-Type M Cassettes. Size range: 4 - 25µm avg. dia. Concentration range: 3,000 - 500,000 cells/ml
For Adherent Cells: Aspirate media and detach cells with appropriate dissociation media (e.g., Accutase).

User Manual:


An electronic copy of the complete Moxi Z user manual is on the USB flash drive that was included in the Moxi Z kit and is available at www.orflo.com.

MoxiChart Software:

An electronic copy of MoxiChart is on the USB flash drive that was included in the Moxi Z kit. MoxiChart facilitates histogram downloading and firmware updates via Bluetooth and is compatible with PC and MAC.

Moxi^Z Mini Automated Cell Counter

1) Turn Moxi ON

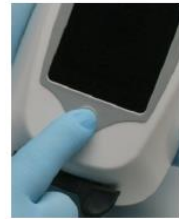2) Depress Tray and Insert Cassette

3) Load 75µL Sample




4) Touch Screen to Start

Touch Here for Cells > 8µm Dia.

Touch Here for Cells < 8µm Dia.

5) Data Summary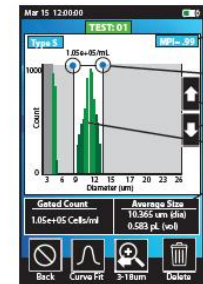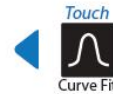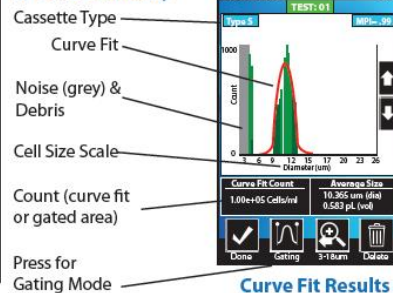


MPI - Moxi Population Index
 Touch to Move Gates
 Adjust Count Scale
 Population Histogram
 Avg. Cell Size (dia. & vol.)
 Press to Change Scale:
 3-18, 3-10,
 3-26µm


Quick Start Guide

Moxi^Z Mini Automated Cell Counter

Saved Histograms: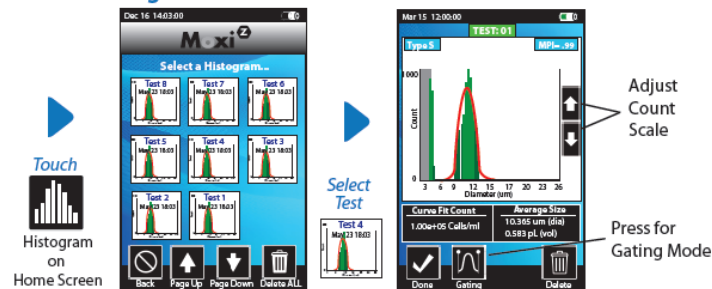

To Obtain Bluetooth ID:

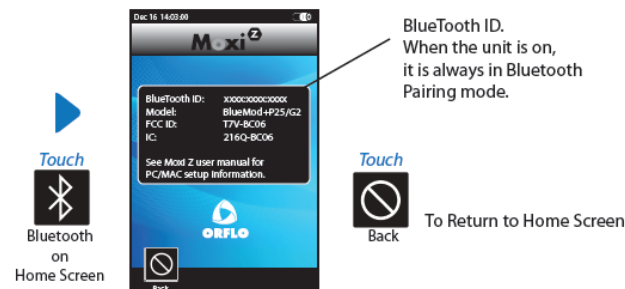To Install MoxiChart Software on the MAC:

- 1) Open ORFLO DISK (on Flash Drive)
- 2) Double Click **MoxiChartInstallMac.dmg**
- 3) Drag **MoxiChart** Icon to Applications folder
- 4) Run MoxiChart by double clicking **MoxiChart** icon.

To Install MoxiChart Software on the PC:

- 1) If prompted to run **MoxiChart Install**, click **YES**. Follow on screen instructions.

OR

- 2) Open ORFLO DISK folder (on Flash Drive) and run **MoxiChart Install**. Follow on screen instructions.